

PHYWE Systeme GmbH & Co. KG
Robert-Bosch-Breite 10
D-37079 Göttingen

Telefon +49 (0) 551 604-0
Fax +49 (0) 551 604-107
E-mail info@phywe.de
Internet www.phywe.com

Operating instructions

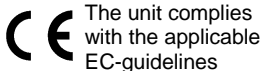


Fig. 1: 12934-01 Cobra SMARTsense EKG

TABLE OF CONTENTS

- 1 SAFETY PRECAUTIONS
- 2 PURPOSE AND CHARACTERISTICS
- 3 FUNCTIONAL AND OPERATING ELEMENTS
- 4 NOTES ON OPERATION
- 5 HANDLING
- 6 TECHNICAL DATA
- 7 SCOPE OF DELIVERY
- 8 ACCESSORIES
- 9 CONFORMITY
- 10 DISPOSAL
- 11 EXPLANATION OF THE SYMBOLS

1 SAFETY PRECAUTIONS



Caution!

- Do not use the Cobra SMARTsense ECG if a device connected to it is also connected to the mains power supply!
- Carefully read these operating instructions completely before operating this instrument. This is necessary to avoid damage to it, as well as for user-safety.
- Only use the instrument for the purpose for which it was designed.
- Only use the instrument in dry rooms in which there is no risk of explosion.
- Protect the instrument from dust, moisture and vapours. Use a slightly moist lint-free cloth to clean the instrument. Do not use aggressive cleaning agents or solvents.
- Take care that no liquid penetrates in through the housing openings, as such penetration would result in damage to Sensor.
- Do not open the unit.

2 PURPOSE AND CHARACTERISTICS

The sensor is used for measuring cardiac activity by way of an ECG (electrophysiology) and for transferring the values to a terminal device, e.g. a tablet computer, smartphone, etc., via Bluetooth or USB.

The sensor is neither suitable nor approved for medical purposes. It is intended solely for didactic purposes. This device must not be used for defined measurements on humans in order to diagnose an illness or disease, i.e. it is not intended

for monitoring, treating or alleviating illnesses or diseases. As a consequence, it is not subject to the strict constraints of the Medical Devices Act.


3 FUNCTIONAL AND OPERATING ELEMENTS

3.1 Operating elements

The sensor has an on-button and two LEDs for indicating the Bluetooth and battery charge status.

On-button 

To switch the sensor on and off in Bluetooth mode, the power button must be pressed for longer than 3s. If the sensor is to be connected via USB, it is not necessary to press the power button.

Bluetooth-LED 

Flashing red every 2 seconds	Not connected
Flashing green every 2 seconds	Connected to the terminal device
Flashing green every 4 seconds	Running measurement

Battery charge LED 

Flashing red every 2 seconds	Low battery
Illuminated red	Active charging process
Illuminated red	Charging process completed

3.2 USB port

The battery, which is permanently installed in the sensor, is charged via the type C USB port. Furthermore, communication with a computer takes place via this interface.

4 NOTES ON OPERATION

For safety reasons, the Cobra SMARTsense sensor may only be used when the sensor is not connected to mains voltage!

To do this, make sure that either

- The Cobra SMARTsense sensor is not connected to a charger via USB.
- If a USB connection to a laptop is established, it is used in battery mode with the power supply disconnected, or
- If there is a USB connection to a standard PC or laptop with a power supply unit, these are connected to the mains voltage via a floating isolating transformer.

This device fulfils all of the technical requirements that are compiled in current EC guidelines. The characteristics of this product qualify it for the CE mark.

This instrument is only to be put into operation under specialist supervision in a controlled electromagnetic environment in research, educational and training facilities (schools, universities, institutes and laboratories).

The individual connecting leads are each not to be longer than 2 m.

The instrument can be so influenced by electrostatic charges and other electromagnetic phenomena (HF, bursts, indirect lightning discharges) that it no longer works within the given specifications. Carry out the following measures to reduce or eliminate the effect of such disturbance: Ensure potential equalization at the PC (especially with Laptops). Use screening. When a total failure of the instrument occurs, unplug it and plug it back in again for a reset.

5 HANDLING

This section describes the start-up of the sensor and the recording of measurement data. Please read this section thoroughly in order to avoid failures or operating errors.

5.1 Charging process

Use a USB-C cable to connect the sensor to a computer or USB charger (not included).

During the charging process, the battery charge LED lights up red. When the charging process is complete, the battery charge LED lights up green. The charging time for a completely discharged battery is 3 hours maximum.



Disconnect the charger at the latest four hours after the completion of the charging process. Otherwise, the service life of the battery may be negatively affected.

5.2 Start-up

Switch the sensor on by pressing the on-button for more than 3 seconds. The Bluetooth LED lights up red. Start the software and select the sensor.

A 9-digit code is printed on the back of the sensor (Fig.2). The last 4 digits of the code are displayed as sensor designation in the software (Fig.3). This enables an exact assignment of the sensors with the software.

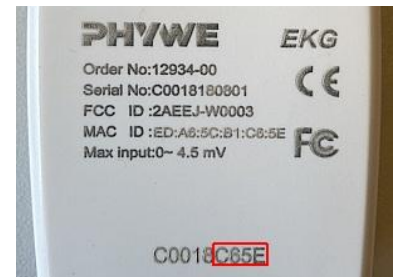


Fig. 2

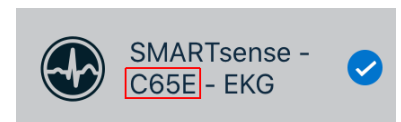


Fig. 3

Selection of the sensor via the Bluetooth interface

Make sure that the Bluetooth interface is activated on the terminal device (PC/Tablet/Smartphone) and that the software is allowed to access the interface.

After the sensor has been selected in the software, the LED flashes green to indicate that the connection has been established correctly. After the sensor has been coupled with the software, the sensor is no longer visible to other users in the software, and therefore can no longer be selected.


If the sensor is switched on and not connected, it switches off automatically after 5 minutes.

Selection of the sensor via the USB interface

For this purpose the sensor must be plugged into the USB port of the end device. It is not necessary to switch on the sensor. The sensor is automatically recognized and displayed. It can be selected and connected directly.

Connecting the electrodes:

Fasten the three disposable electrodes in place: on the inside of both wrist and on the inside of the left lower leg, right above the ankle. Connect the three colour-coded individual measuring leads as follows:

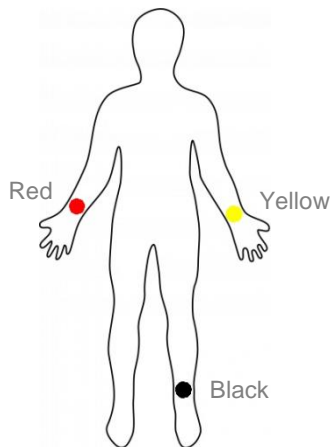

Then, plug the jack plug into the front socket of the sensor.

5.3 Recording of measurement data

During the measurement, it should be taken into account that even slight muscle activity can influence the measurement. This is why the measurement must be performed while sitting or lying down in a relaxed manner. The person should place their hands on their thighs during the measurement.

Adjust a data rate of 200 Hz minimum in the software and start the measurement.

6 TECHNICAL DATA

Operating temperature range: 5 - 40°C

Rel. humidity < 80%

Measuring range	0...4.5 mV
Resolution	4.5 µV

Max. data rate	1000 Hz
Battery capacity	250 mAh
Max. wireless range (open field)	30 m
Dimensions (length x width x height)	80 x 40 x 23 mm
Weight	40 g

7 SCOPE OF DELIVERY

The extent of delivery is as follows

- Cobra SMARTsense EKG 12934-01
- USB cable type C 07935-00
- 6 EKG electrodes and connecting cable
- Operating instructions

8 ACCESSORIES

The following accessories are available:

- Cobra SMARTlink 12999-99
- USB-charger 07934-99
- USB cable type C 07935-00
- Set with 20 Selfsticking Electrodes 12929-00
- USB-Bluetooth-Adapter 07936-00
- Software measureLAB 14580-61
- Free measureApp available from supplier portals

iOS



Android



Windows



9 CONFORMITY



PHYWE Systeme GmbH & Co.KG hereby declares that the radio system type 12934-01 complies with the 2014/53/EU directive. The complete text of the EC Declaration of Conformity is available at the following Internet address:

www.phywe.com/en/ec-declaration

10 DISPOSAL

The packaging mainly consists of environmentally-friendly materials that should be returned to the local recycling stations.



Do not dispose of this product with normal household waste. If this unit needs to be disposed of, please return it to the address that is stated below for proper disposal

PHYWE Systeme GmbH & Co. KG
Abteilung Kundendienst
Robert-Bosch-Breite 10
D-37079 Göttingen

Telefon	+49 (0) 551 604-274
Fax	+49 (0) 551 604-246

11 EXPLANATION OF THE SYMBOLS



Attention!
Potentially damaging situation (damage to property)
General danger area
Observe operating instructions



Note
Important information about the device
Observe operating instructions