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GENERAL DESCRIPTION

All the appliance is made in metal:

- The burner, which, depending on the nozzle, is suitable for operation with all types of gas (Town Gas - L.P.G. Gas - Natural Gas, is made of chromate brass.
- The Bunsen Burners are equipped with Cock with Safety Needle provided with an anti-withdrawing locking device.
- "SEKUR FLAM" and "SEKUR FLAM POLIGAS" Bunsen Burners are provided with Thermocouple-controlled and Thermoelectric Safety Valve to cut OFF the gas supply in case of accidental extinguishment of the flame.
- The valves have been designed, built and controlled according to the essential safety requirements of directive 2014/68 / UE (PED).
- The new line Bunsen will be made with a steel plated disk, painted by an epoxy and Acidproof painting. We also added a PVC ring to not scratch the surfaces.

START THE OPERATION

Do not start to use the product before reading directions for use.

Warn of Danger

To prevent any accidents caused by improper use of the product, please follow the following safety rules

SAFETY RULES

The installation of the burner to the gas should preferably be performed by a licensed contractor.

Before installing the burner to the gas network to ensure that the gas pressure get it right.

Make sure before you connect the burner to the network that the tap screw is closed.

Make sure the inlet is properly connected to the hose and that there are no gas leaks.

Do not connect the burner to gas, without mounting a pressure reducer.

Do not connect to the Bunsen gas directly to gas container.

Do not change any part of the Bunsen (holes, passages, nozzles etc.)..

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SPECIFICATIONS

All types of Bunsen burner built by us (Easy-Meker-Teclu-Poligas):

- ***SINGLE FAMILY TYPE***
 - Easy - operated by: Natural Gas, Liquid Gas, Natural Gas.
 - Easy - Poligas: works with all types of Gas.
- ***FAMILY TYPE “SEKUR FLAM”***
 - Sekur-Flam - working with: Natural Gas, Liquid Gas, Natural Gas.
- ***FAMILY TYPE “SEKUR FLAM-POLIGAS”***
 - Poligas - working with all types of Gas.

They are supplied with a sheet that reports the technical specifications and certifies their testing.

HOW TO USE

BUNSEN BURNER ON SIMPLE

- Turn the compass air entry closing completely;
- Turn in one turn the knob on the gas inlet valve;
- There is placed on top of the Bunsen burner, with an open flame or with a piezoelectric lighter, causing the ignition;
- Proceed with gradual opening up of gas and air until they reach desired flame.

BUNSEN BURNER IGNITION “SEKURFLAM”

- Turn the compass air entry closing completely;
- Turn in one turn the knob on the gas inlet valve, thermoelectric valve does not allow the entry of gas;
- Press spring valve will exit from the Gas Superior of the Bunsen burner;
- There is placed on top of the Bunsen burner, with an open flame or with a piezoelectric lighter, causing the ignition;
- Hold the valve button for 10 seconds to allow the thermocouple to reach operating temperature fully open.

INTERVENTION OF THERMOELECTRIC SAFETY VALVE

If you were to accidentally turn off the heat, in a time of about 40 seconds due to the lowering of the temperature of the thermocouple spout located near the flame, the thermoelectric valve, through the built-in magnetic core, automatically closes the gas outlet.

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ON BUNSEN “SEKURFLAM” POLIGAS

- *These lamps are equipped with special adjustable nozzle that can be used with any type of gas;*
- *Turn the compass dial entry or air to close completely;*
- *Turn in one turn the knob on the faucet or the bell of the burner tube for gas inlet, the valve does not allow the entrance of the thermal gas;*
- *Press the button on the valve spring causes the release of gases from the Superior of the Bunsen burner;*
- *There is placed on top of the Bunsen burner, with an open flame or with a piezoelectric lighter, causing the ignition;*
- *Hold the valve button for 10 seconds to allow the thermocouple to reach operating temperature fully open;*
- *Proceed with gradual opening of the gas through the knob, pipe, or the bell. REGULAR AIR, with the compass or the ring under the bell until you reach the desired flame.*

INTERVENTION OF THERMOELECTRIC SAFETY VALVE

If you were to accidentally turn off the heat, in a time of about 40 seconds sag due to the temperature of the tip of the thermocouple in the vicinity of the flame, the thermal valve through the magnetic core incorporated automatically closes the gas outlet.